



**Thermo**  
ELECTRON CORPORATION  
*Karl Fischer Titrator*

Log # **KF1086**

**Overview**

---

**Sample** POLYACRYLAMIDE GRANULES

**Industry** POLYMERS AND PLASTICS

**Group Number** 27

**Typical moisture** 7%

**Sample size** 0.7g

**Summary of  
procedure**

Method Parameters

<b>Instrument</b>	Turbo 2 Volumetric Karl Fische	<b>Titration Current</b>	
<b>Volumetric Reagent</b>	Reidel de Haen Composite 5	<b>Endpt Rate</b>	
<b>Volumetric Solvent</b>	MEOH CONH2 1/1	<b>Stirrer/blender speed</b>	3
<b>Pump Tubing</b>	SILICONE	<b>Extract time</b>	15.0min
<b>Vessel/Vessel Tubing</b>	STD 750mL	<b>Mode of analysis</b>	SAMPLE ONLY
<b>Endpt/Step Level</b>	8/5	<b>Vaporizer oven</b>	
<b>Endpt Time</b>	10sec	<b>Result units</b>	% wt/wt
<b>Anode Reagent</b>		<b>Source and date</b>	TW (075)
<b>Cathode Reagent</b>		<b>Cross-reference log</b>	KF1035

Results

<b># of Trials</b>	3	<b>Mean Value</b>	6.95%	<b>C. V.</b>	0.7%
--------------------	---	-------------------	-------	--------------	------

**Comments**