

Application Notes for Protein in Milk-Whey-Soy Powders


Summary

Product: milk-whey-soy
Description: off-white, fine powder
Parameter Measured: protein
Typical Range: 10-25%
Typical RMS Error: 0.2-0.6%

Application Notes:

General: This application note covers a variety of powder types and formulations.
Presentation: Off-line over a turntable, on-line over a transport conveyor.
Benefits: Protein content needs to be tightly controlled.

Data Plot:

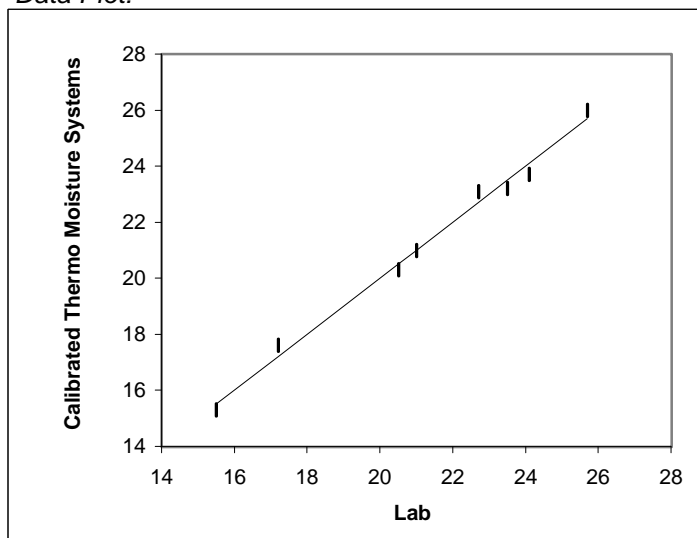

Cal. Data:

Sample	% Protein	
	Lab	TMS
1	15.50	15.3
2	17.20	17.6
3	20.50	20.3
4	21.00	21.0
5	22.70	23.1
6	23.50	23.2
7	24.10	23.7
8	25.70	26.0