

DYNAMIC CONTACT ANGLE ACCESSORY FOR THIN FILM SUBSTRATES

The ability to measure the wetting properties of thin films by the Wilhelmy Plate Technique has been a challenge due to the lack of rigidity in many film substrates.

Polymer film substrates such as Mylar are being used today in many industrial applications including many packaging and printing operations. These materials are typically prepared as very thin film sheets with an average thickness of less than 15 microns. The surfaces of these substrate materials are often subjected to dip coating operations to control the wettability of the surface to fit the desired end-use application.

Early attempts to characterize the wettability of these thin film substrates by the Wilhelmy Plate Immersion Technique were unsuccessful as the film sample (typically 5-20 microns thick, cut 10-30 millimeters square and held in place by a traditional single point contact sample holder) passed through the air/liquid interface, the sample was observed to lay down or bend, failing to penetrate the interface perpendicular to the liquid surface layer (see *Figure 1*). This phenomenon was first observed with a 12 micron thick sample of Mylar film prepared in the manner described above.

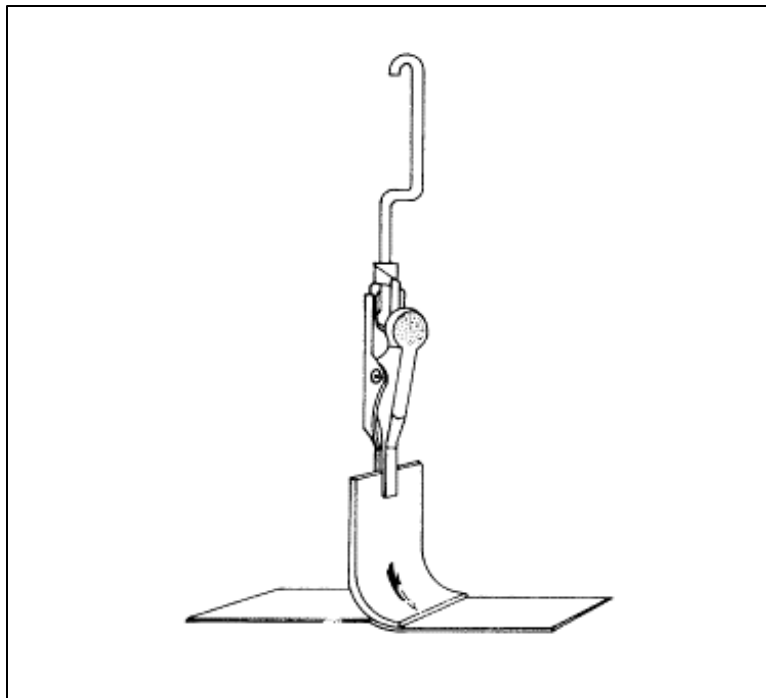


Figure 1: Mylar film held by traditional single point sample holder fails to penetrate the air/liquid interface.

After studying this phenomenon and experimenting with several options aimed at overcoming the flotation, Thermo Cahn application engineers came up with a solution. A new type of sample holder was developed (see *Figure 2*) to increase the rigidity of the film by forcing the sample into a slight curved orientation perpendicular to the air/liquid interface. The Thermo Cahn Thin Film Sample Holder (*Part #21107-02*) is recommended for use with most thin non-rigid materials including a variety of thin films and foils. Samples are prepared in the usual manner with vertical sides approximately 10 millimeters apart (sample holder 12 millimeters wide) and inserted into the thin film sample holder with the length of sample exposed for wettability scanning limited to 5-10 mm depending on the flexibility of the sample.

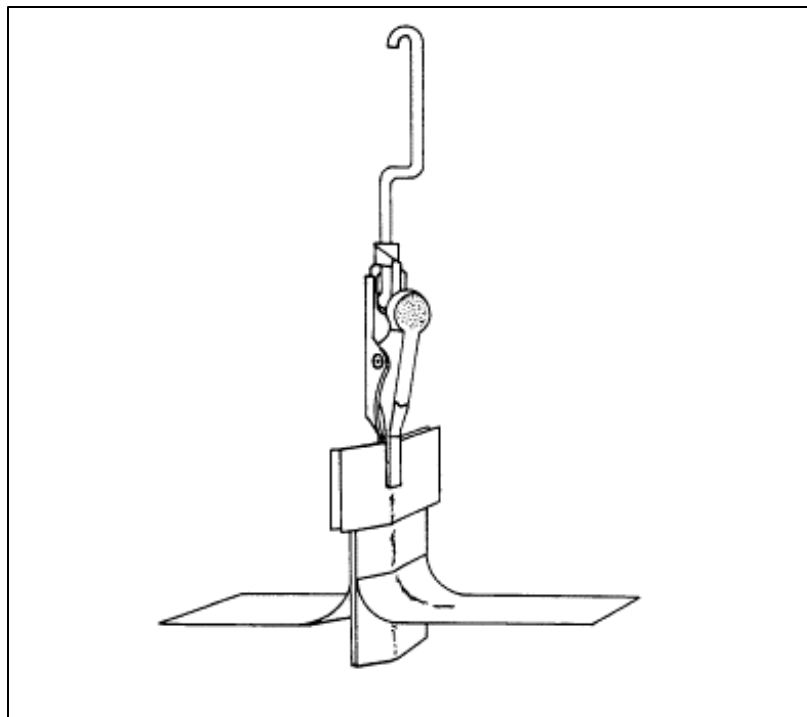


Figure 2: Mylar film readily penetrates air/liquid interface when held rigidly between copper-beryllium plates of Cahn Thin Film

The holder is constructed of a rigid alligator clip design with two parallel copper-beryllium plates attached at one end. The plates have matched curvatures to sandwich the film sample uniformly and thus force the film into a rigid conformation. All users of Thermo Cahn DCA Systems can utilize the Thin Film Sample Holder.

Thermo Cahn's Thin Film Sample Holder opens up a whole new class of materials to experimentation by the Dynamic Wilhelmy Plate Method.

Thermo Material Characterization

A Thermo Electron Company

International
Thermo Haake
Dieselstr. 4 · 76227 Karlsruhe
Germany
Tel. +49(0)721 40 94-0
Fax +49(0)721 40 94-300
Hotline +49(0)18050 HAAKE
info@thermohaake.com
www.thermohaake.com

USA
Thermo Haake
5225 Verona Road
Madison, WI 53711
Tel. 608-327-6777
Fax 608-273-6827
infousa@thermohaake.com
www.thermohaake.com

France
Thermo Rheo
99 route de Versailles
91160 Champlan
Tel. +33(0)1 64 54 01 01
Fax +33(0)1 64 54 01 87
info@thermorheo.com
www.thermorheo.com

UK
Thermo PRISM
Emerald Way
Stone Business Park, Stone
Staffordshire ST15 0SR
Tel.+44(0)1785 813 648
Fax+44(0)1785 814 633
info@thermoprism.com
www.thermoprism.com