

Thermo Scientific Orion 15 mm Electrode Holder for the Swing Arm Electrode Stand

General Information

The Thermo Scientific Orion swing arm electrode stand, Cat. No. 090043, is assembled and shipped with an electrode holder that is designed to hold up to five electrodes with 12 mm body diameters. See **Figure 1**.

An additional electrode holder, Cat. No. 090044, can be purchased separately and is designed to hold an electrode with a 15 mm body diameter. This electrode holder can fit one 15 mm diameter electrode and up to four 12 mm diameter electrodes. See **Figure 2**. The replacement electrode holder gently grips the 15 mm (0.60 inch) diameter electrode body, which keeps the electrode stable in a beaker or other sample container.

Figure 1
12 mm Electrode Holder
Supplied with the Stand


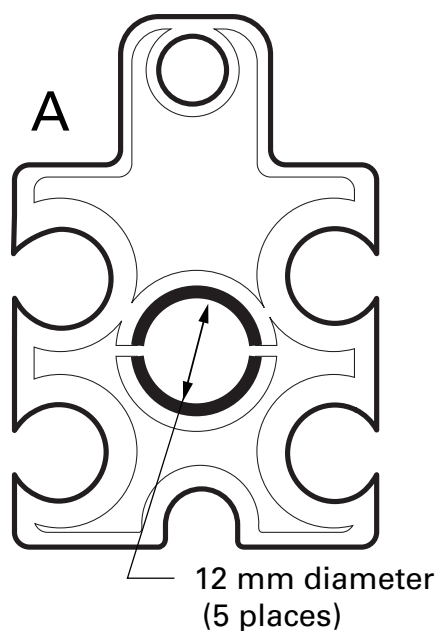



Figure 2
15 mm Electrode Holder

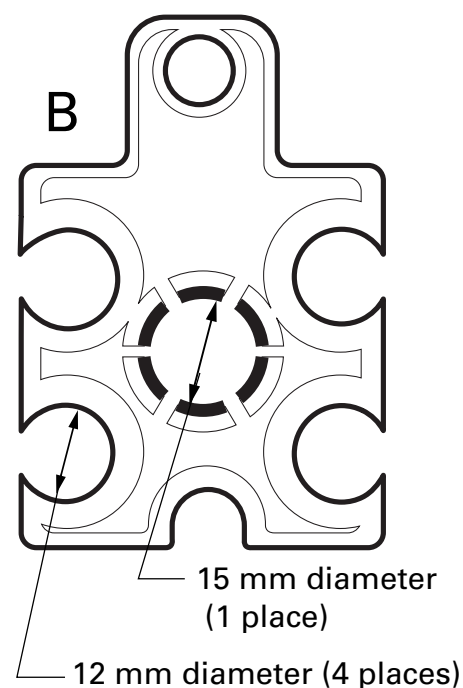

Assembly Instructions

1. Remove the pin that connects the swing arm and electrode holder.
2. Remove the existing electrode holder.
3. Install the 15 mm electrode holder.
4. Replace the pin that connects the swing arm and electrode holder.

Figure 3
15 mm Electrode Holder
Installation on Stand

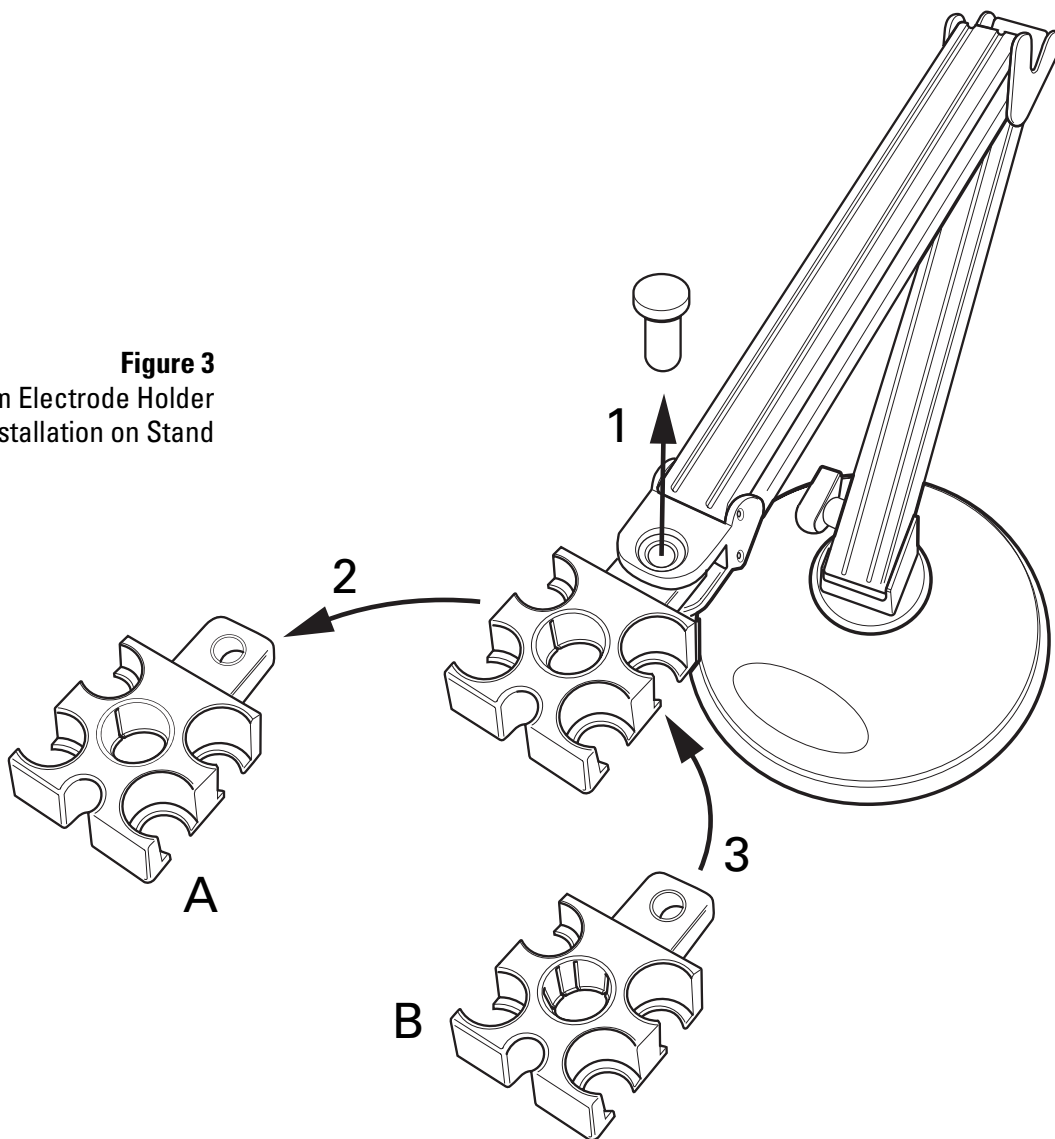

Ordering Information

Cat. No.	Description
090043	Swing arm electrode stand, includes 12 mm electrode holder
090044	15 mm electrode holder for use with swing arm electrode stand

©2009 Thermo Fisher Scientific Inc. All rights reserved. All trademarks are the property of Thermo Fisher Scientific Inc. and its subsidiaries.



Water Analysis Instruments

North America
166 Cummings Center
Beverly, MA 01915 USA
Toll Free: 1-800-225-1480
Tel: 1-978-232-6000
Dom. Fax: 1-978-232-6015
Int'l Fax: 978-232-6031

Europe
P.O. Box 254, 3860 AG Nijkerk
Wallerstraat 125K, 3862 CN
Nijkerk, Netherlands
Tel: (31) 033-2463887
Fax: (31) 033-2460832

Asia Pacific
Blk 55, Ayer Rajah Crescent
#04-16/24, Singapore 139949
Tel: 65-6778-6876
Fax: 65-6773-0836

www.thermo.com/water

255024-001 Rev. A 05-09

Thermo
SCIENTIFIC