






H-6000A
 Stainless Steel Rotor Body
 Aluminum Buckets
 6-Place, 1000 ml
 Tube Dimensions:
 98 × 175 mm
 Swinging-Bucket
 Maximum Speed:
 5,000 rpm
 Rotor Mass: 24.95 kg

Buckets	RCF (g)	R (cm)	Speed (rpm)
Max.	7,277	26.06	5,000
Microplate Carriers	RCF (g)	R (cm)	Speed (rpm)
Max.	5,186	18.57	5,000

For use in **RC3BP®**, **RC-3C Plus**, **RC-3C**,¹ **RC-3B Plus** and **RC-3B**¹ centrifuges

Catalog No. 11250	Catalog No. 11320	Accessories
Includes: H-6000A Rotor Rotor Cover 11237 Rotor Buckets, Anodized Aluminum, (3 pkgs, 2/pkg) 44040 Rotor Locking Tool 11246 Rotor Cleaning Kit 12259 Balancing Discs, Rubber (1 pkg, 6/pkg) 00335 Instruction Manual 11243	Includes: H-6000A Rotor includes all accessories in catalog no. 11250 <u>without buckets</u>	Microplate Carriers (set of 2) 11267 Rotor Extractor 11445 Sealing Covers (set of 2) High Heat Polycarbonate for threaded buckets 11776 Adapter Removal Post 00392 Sealing Cover O-rings (set of 6) 12995 Dust cover nonsealing O-ring 66721 (1)
 00335		 11776
		 11267

TUBES, BOTTLES

NOTE: HBB-6 Conversion Kit **11758** (includes (6) HBB-6 buckets, manual, disks)

CAPACITY (ml)		DESCRIPTION	MAX SPEED (rpm)	DIMENSIONS (mm)	TUBE CAT. NO. (Qty/Pkg)	SEALING ASSEMBLY*		ADAPTER REQ'D		NOTES
NOMINAL	FILL					CAT. NO. (Qty/Pkg)	DESCRIPTION	CAT. NO. (Qty/Pkg)	NO. OF PLACES	
1,000	960	Polycarbonate Oak Ridge Bottle	5,000	98 × 175	03947 (4)	03280 (12)	Polypropylene, Sealing	—	—	Nonsealing Cap 03928 (12/pkg)
1,000	1,000	Polypropylene Bottle	5,000	98 × 175	03695 (4)	Included with Bottle	Polypropylene, Nonsealing	—	—	Replacement Nonsealing Cap 00707 available
500	—	Polypropylene Corning Tissue Culture Bottle, Conical	—	96 × 146	431123 ²	Included with Bottle	Polypropylene, Plug	431124 ²	1	
450 to 550	—	Blood Bag, Triple, Quad	—	—	—	—	—	00512 (1)	1	
450	—	Blood Bag, Single and Double	—	—	—	—	—	00511 (1)	1	
500	425	Polycarbonate Oak Ridge Bottle	5,000	70 × 160	03944 (6)	03280 (12)	Polypropylene, Sealing	00444 (1) 00511 (1)	1 1	Adapters 00444 and 00511 are both required. Nonsealing Cap 03928 (12/pkg) available.
500	425	Polypropylene Oak Ridge Bottle	5,000	70 × 160	03943 (6)	03280 (12)	Polypropylene, Sealing	00444 (1) 00511 (1)	1 1	Adapters 00444 and 00511 are both required. Nonsealing Cap 03928 (12/pkg) available.
315	360	Stainless Steel Bottle	5,000	61 × 153	00522 (1)	00518 (1)	Stainless Steel, Sealing	01092 (1)	1	Requires Wrench 01014 to seal bottle.
285	330	Stainless Steel Bottle	5,000	61 × 140	00530 (1)	Included with Bottle	Stainless Steel, Sealing	001092 (1)	1	

⁽¹⁾Instrument no longer available.

⁽²⁾Available from Corning CoStar. In U.S. call 800-492-1110.

*Sealing Assembly replacement parts may be available. See section E for availability.

(continued on next page)

TUBES, BOTTLES FOR H-6000A (continued)

CAPACITY (ml)		DESCRIPTION	MAX SPEED (rpm)	DIMENSIONS (mm)	TUBE CAT. NO. (Qty/Pkg)	SEALING ASSEMBLY*		ADAPTER REQ'D		NOTES
NOMINAL	FILL					CAT. NO. (Qty/Pkg)	DESCRIPTION	CAT. NO. (Qty/Pkg)	NO. OF PLACES	
250	250	Polypropylene Oak Ridge Bottle	5,000	61 × 124	03937 (12)	03278 (12)	Polypropylene, Sealing	01092 (1)	1	Nonsealing Cap 03927 (12/pkg) available.
250	265	Polycarbonate Oak Ridge Bottle	5,000	61 × 124	03939 (12)	03278 (12)	Polypropylene, Sealing	01092 (1)	1	Nonsealing Cap 03927 (12/pkg) available.
250	—	Tissue Culture Bottle, Conical	—	60 × 177	—	—	—	01092 (1) 03095 (1)	1	Both adapters are required.
200	200	Nunc Tissue Culture Bottle, Conical	4,900	60 × 135	03094 (4)	Included with Bottle	Polypropylene	01092 (1) 03095 (1)	1 1	Both adapters are required.
200	350	Stainless Steel Tube	5,000	61 × 124	00521 (1)	—	—	01092 (1)	1	
150	170	Corex® Bottle (No longer available)	5,000	53 × 113	03229 (1)	Included with Bottle	Polypropylene, Nonsealing	00372 (1) 01092 (1)	1 1	Both adapters are required.
125	150	Corex® Bottle (No longer available)	5,000	53 × 118	03231 (1)	Included with Bottle	Polypropylene, Nonsealing	00372 (1) 01092 (1)	1 1	Both adapters are required.
100	134	Pyrex® Bottle	3,000	44 × 137	03230 (1)	Included with Bottle	Polypropylene, Nonsealing	00371 (1) 01092 (1)	1 1	Both adapters are required.,
50	Accepts all Tubes, Bottles, and Adapters for SS-34 and SA-600 Rotors; requires Adapter 00894 (1/pkg, 5-place) PLUS any adapter specified in SS-34/SA-600 tube chart. See page pages B-30 and B-34.									
50	—	Tissue Culture Tube, Conical	—	30 × 115	—	—	—	00436 (1) or 00445 (1)	5 or 3	Both adapters accommodate freestanding and conical tissue culture tubes.
35	—	Polyallomer Nalgene® Conical Oak Ridge Tube	—	29 × 105	—	—	—	00436 (1) or 00445 (1)	5 or 3	Both adapters accommodate freestanding and conical tissue culture tubes.
35	—	Polycarbonate Nalgene® Conical Oak Ridge Tube	—	29 × 105	—	—	—	00436 (1) or 00445 (1)	5 or 3	Both adapters accommodate freestanding and conical tissue culture tubes.
16	16	Polypropylene Flanged Tube	5,000	18 × 100	03244 (50)	03299 (50)	Polypropylene, Snap-on	00892 (1)	12	
16	17	Polycarbonate Flanged Tube	5,000	18 × 100	03243 (50)	03269 (50)	Polypropylene, Snap-on	00892 (1)	12	
15	—	Urinalysis Tube, Conical	—	17 × 120	—	—	—	00892 (1)	12	
15	—	Tissue Culture Tube, Conical	—	17 × 120	—	—	—	00447 (1)	8	
15	—	Vacutainer® Tube	—	16 × 127	—	—	—	00385 (1)	18	
12	14	Pyrex® Tube	4,000	18 × 102	03109	—	—	00894 (1)	5	5 Rubber Adapters 00363 required with 00894 .
15	18	Corex® Tube	4,000	18 × 102	03286 (25)	—	—	00892 (1)	12	Rubber Pads 03667 (12/pkg) recommended.
12	12	Polypropylene Flanged Tube	5,000	16 × 100	03116 (50)	03266 (50)	Polypropylene, Snap-on	00385 (1)	18	
12	12	Polycarbonate Flanged Tube	5,000	16 × 100	03115 (50)	03266 (50)	Polypropylene, Snap-on	00385 (1)	18	

*Sealing Assembly replacement parts may be available. See section E for availability.

(continued on next page)

All trademarks are the property of their respective owners.

TO ORDER, CALL 1-800-522-SPIN (7746)

TUBES, BOTTLES FOR H-6000A (continued)

CAPACITY (ml)		DESCRIPTION	MAX SPEED (rpm)	DIMENSIONS (mm)	TUBE CAT. NO. (Qty/Pkg)	SEALING ASSEMBLY*		ADAPTER REQ'D		NOTES
NOMINAL	FILL					CAT. NO. (Qty/Pkg)	DESCRIPTION	CAT. NO. (Qty/Pkg)	NO. OF PLACES	
10	—	Vacutainer® Tube	—	16 × 100	—	—	—	00385 (1)	18	
10	11	Polyallomer Thin-Walled Flanged Tube	5,000	13 × 100	03124 (50)	—	—	00384 (1)	22	
10	—	Glass Tube	—	13 × 100	—	—	—	00384 (1)	22	
10	—	Pyrex® Tube	—	13 × 100	—	—	—	00384 (1)	22	
10	9.1	Polypropylene Oak Ridge Bottle	4,000	16 × 83	03929 (50)	03279 (25)	Polypropylene, Sealing	00894 (1) and 00402 (5)	5	Nonsealing Cap 03924 (25/pkg) available.
10	9.1	Polycarbonate Oak Ridge Bottle	4,000	16 × 83	03020 (25)	03279 (25)	Polypropylene, Sealing	00894 (1) and 00402 (5)	5	Nonsealing Cap 03924 (25/pkg) available.
7	—	Vacutainer® Tube	—	13 × 100	—	—	—	00384 (1)	22	
7	7	Polypropylene Flanged Tube	5,000	13 × 100	03121 (50)	03265 (50)	Polypropylene, Snap-on	00384 (1)	22	
7	7	Polycarbonate Flanged Tube	5,000	13 × 100	03120 (50)	03265 (50)	Polypropylene, Snap-on	00384 (1)	22	
5	6	Pyrex® Tube	5,000	12 × 75	03102 (50)	—	—	00857 (1) 00858 (1)	33 33	Top Level Adapter 00858 must be used with Bottom Level Adapter 00857 . Adapter 00857 may be used alone.
5	—	Standard Glass Tube	—	12 × 75	—	—	—	00857 (1) 00858 (1)	33 33	Top Level Adapter 00858 must be used with Bottom Level Adapter 00857 . Adapter 00857 may be used alone.
5	—	Vacutainer® Tube	—	10.25 × 103	—	—	—	00857 (1)	33	
5	—	Vacutainer® Tube	—	13 × 75	—	—	—	00857 (1)	33	Only 14 tubes per adapter due to size of plug.
4	—	Vacutainer® Tube	—	10.25 × 82	—	—	—	00857 (1)	33	
4	—	Standard Glass Tube	—	10 × 75	—	—	—	00857 (1) 00858 (1)	33 33	Top Level Adapter 00858 must be used with Bottom Level Adapter 00857 . Adapter 00857 may be used alone.
4	4	Polypropylene Flanged Tube	5,000	11 × 75	03105 (50)	03264 (50)	Polypropylene, Snap-on	00857 (1) 00858 (1)	33 33	Top level Adapter 00858 must be used with Bottom Level Adapter 00857 . Adapter 00857 may be used alone.
4	4	Polycarbonate Flanged Tube	5,000	11 × 75	03104 (50)	03264 (50)	Polypropylene, Snap-on	00857 (1) 00858 (1)	33 33	Top level Adapter 00858 must be used with Bottom Level Adapter 00857 . Adapter 00857 may be used alone.
3	3	Pyrex® Tube	5,000	10 × 75	03100 (50)	—	—	00857 (1) 00858 (1)	33 33	Top Level Adapter 00858 must be used with Bottom Level Adapter 00857 . Adapter 00857 may be used alone.
1.5	—	Microtube	—	—	—	—	—	04871 (1) 00857 (1)	— 33	Adapter plate 04871 must be used together with adapter 00857 .
—	—	Microplate, 3 Standard or 1 Deepwell	—	—	—	—	—	—	3 1	Carrier 11267 is used in place of bucket.

*Sealing Assembly replacement parts may be available. See section E for availability.

All trademarks are the property of their respective owners.