

Thermo Scientific Orion Star Series Printer

The Star Series inkjet printer ships with one black ink cartridge only.

Thermo Scientific power cords for US, UK, Europe and Australia have been included with the printer. The AC adapter is universal and will work with most AC voltages.

Before operating the printer, DIP switch SW1-3 must be set to the OFF position and DIP switch SW1-4 must be set to the ON position. Setting DIP switch SW1-3 to the OFF position disables the paper cutter when a self-test is performed on the printer. Setting DIP switch SW1-4 to the ON position instructs the printer that there is only one ink cartridge installed.

Caution: If the printer is not turned off when the DIP switch cover is removed, an electrical short may result that can damage the printer.

To change the DIP switch settings

1. Turn the printer off.
2. Remove the screw from the DIP switch cover and then remove the DIP switch cover from the printer. The DIP switch cover is located on the bottom of the printer.
3. Set the DIP switches using a pointed tool, such as tweezers or a small screwdriver.
4. Replace the DIP switch cover and then secure it with the screw.
5. Turn the printer on.

Ordering Information

Cat. No.	Description
1010006	Star Series inkjet printer with cable
1010008	Ink cartridge for the Star Series printer
1010009	Paper for the Star Series printer

Environmental Instruments Water Analysis Instruments

166 Cummings Center
Beverly, MA 01915 USA
Toll Free: 1-800-225-1480
Tel: 1-978-232-6000
Dom. Fax: 1-978-232-6015
Int'l Fax: 978-232-6031

www.thermo.com/water

© 2007 Thermo Fisher Scientific Inc.
All rights reserved.