

Lithium isotope measurements using Finnigan™ NEPTUNE Multicollector ICP-MS

Claudia Bouman, Johannes B. Schwieters, Thermo Electron (Bremen) GmbH, Bremen, Germany

Key Words

- Finnigan™ NEPTUNE
- International Rock Standards
- Lithium Isotopes
- Multicollector ICP-MS
- NIST L-SVEC
- Static Collection

Introduction

Precise isotopic analysis of Li is required in a number of different applications, including geochemistry and nuclear sciences. Thermal ionization mass spectrometry (TIMS) is the preferred technique used for precise measurements of the ${}^7\text{Li}/{}^6\text{Li}$ ratio. A new technique for Li isotope analysis is Multicollector Magnetic Sector Inductively Coupled Plasma Mass Spectrometry (MC-ICP-MS) (e.g. Tomascak et al. 1999a,b). Here, we report on the short-term reproducibility of Li isotope analyses of NIST L-SVEC lithium carbonate standard and two international rock standards, BHVO-1 (USGS) and JG-2 (GSJ) with the Finnigan NEPTUNE MC-ICP-MS.

Experimental

Samples are introduced in the ICP using a low flow self-aspirating PFA concentric nebulizer (uptake ~ 80 $\mu\text{l}/\text{min}$) and a quartz dual spray chamber ("stable introduction system"). The Finnigan NEPTUNE MC-ICP-MS from Thermo Electron Corporation, allows simultaneous measurement of ${}^6\text{Li}^+$ and ${}^7\text{Li}^+$ using the moveable low mass and high mass Faraday cups of the collector array (on platforms L4 and H4, respectively). The analysis sequence was: blank, L-SVEC standard, blank, sample 1, blank, L-SVEC standard, blank, sample 2, blank etc., so that a blank (an 'on peak' baseline) was measured before and after each sample or standard. 0.5% Triton X100 (Merck)

was added to standards, samples and blanks to improve beam stability. After initial uptake of sample (60 s), the analysis was started. The reproducibility of Li isotope measurements was evaluated by analyzing a 200 ppb Li L-SVEC standard giving a typical ${}^7\text{Li}$ beam intensity of ~4 V (4×10^{11} A at 10^{11} Ω resistor). With an uptake rate of 80 $\mu\text{l}/\text{min}$, approximately 80 ng of sample were consumed during one analysis lasting about 5 minutes (3 blocks of 10 ratios, 8 s per ratio measurement). To minimize memory effects, the aspiration system was rinsed for at least 5 minutes in 3% HNO_3 after each analysis to ensure that the background reached an appropriate baseline value.

Results

L-SVEC Li standard

Figure 1 shows the short-term reproducibility of the Li isotope measurements on the L-SVEC standard solution. One analysis of three blocks (30 ratios) results in an internal precision of ~ 0.2 ‰ (1 σ). The external reproducibility (n = 8) for ${}^7\text{Li}/{}^6\text{Li}$ is also about 0.2 ‰ (1 σ). While the mass bias is large, around 25 % per amu (${}^7\text{Li}/{}^6\text{Li}$ L-SVEC = 12.02 ± 0.03 ‰; Flesch et al., 1973 vs. 15.0-15.2 ‰, Figures 1 and 2) and varies with time, long-term variations are rather small (within 2 ‰) and the ${}^7\text{Li}/{}^6\text{Li}$ ratio of the L-SVEC standard varies around a mean of 15.

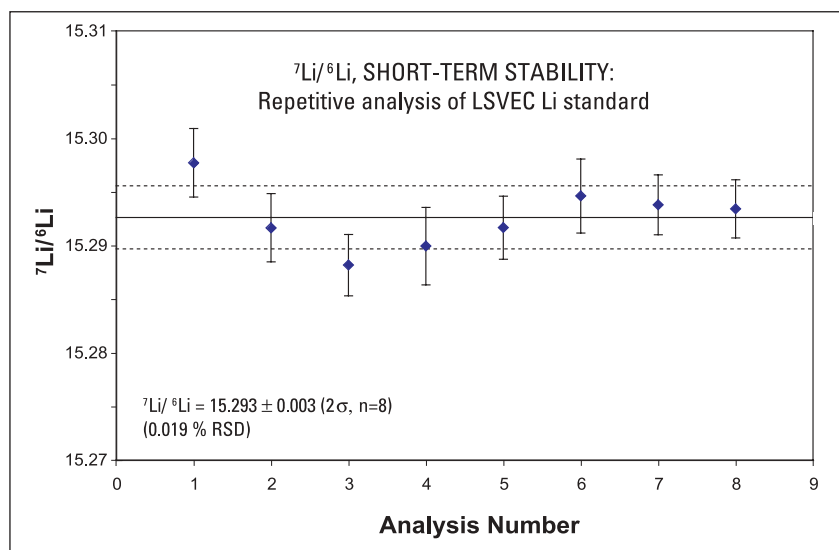


Figure 1: Short-term reproducibility of Li isotope measurements (L-SVEC Li standard, 200 ppb). Duration: 4 hours.

International rock standards

Li isotope analyses of two international standards, BHVO-1 and JG-2, were performed to ensure the reliability of Li isotope data obtained by MC-ICP-MS. The measured ${}^7\text{Li}/{}^6\text{Li}$ of the samples were normalized to the mean of the measured ${}^7\text{Li}/{}^6\text{Li}$ of the standard runs before and after (linear interpolation). Data are presented as deviations of the measured ${}^7\text{Li}/{}^6\text{Li}$ in part per thousand from the ${}^7\text{Li}/{}^6\text{Li}$ of the standard ($\delta^7\text{Li}$).

$$\delta^7\text{Li} (\text{‰}) = \left(\frac{R_{\text{sa}}}{R_{\text{st}}} - 1 \right) \times 10^3$$

where R_{sa} is the ${}^7\text{Li}/{}^6\text{Li}$ ratio of the sample and R_{st} is the ${}^7\text{Li}/{}^6\text{Li}$ ratio of the L-SVEC standard.

Figure 2 shows the results of the Li isotope analyses of BHVO-1 and JG-2. The Li isotope compositions for JG-2 are in good agreement with the values reported by Oi et al. (1997) and James and Palmer (2000). There are no published data for BHVO-1. However, our results for BHVO-1 agree well with MC-ICP-MS analysis for a suite of other Hawaiian basalts, which define a range in $\delta^7\text{Li}$ of +3.0 to +4.8‰ (Tomascak et al., 1999b).

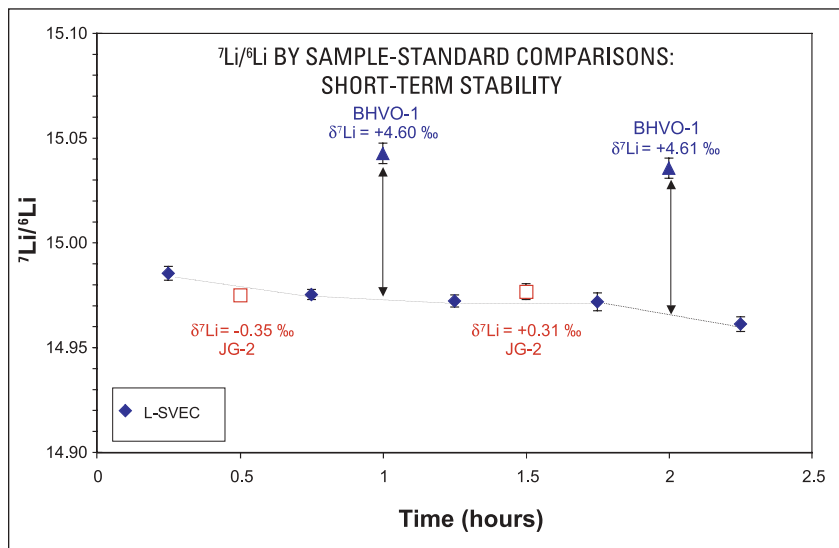


Figure 2: Li isotope analyses of international rock standards BHVO-1 and JG-2. δ values (in ‰) are given relative to the L-SVEC Li standard.

References

- Flesch G., Anderson A.R., and Svec H.J. (1973) A secondary isotopic standard for ${}^6\text{Li}/{}^7\text{Li}$ determinations, *International Journal of Mass Spectrometry and Ion Physics* 12 265-272.
- James R.H., and Palmer M.R. (2000) The lithium isotope composition of international rock standards, *Chemical Geology* 166 319-326.
- Oi T., Odagiri T., and Nomura M. (1997) Extraction of lithium from GSJ rock reference samples and determination of their lithium isotopic composition, *Analytica Chimica Acta* 340 221-225.
- Tomascak P., Carlson R.W., and Shirey S.B. (1999a) Accurate and precise determination of Li isotopic compositions by multi-collector sector ICP-MS, *Chemical Geology* 158 145-154.
- Tomascak P., Tera F., Helz R.T., and Walker R.J. (1999b) The absence of lithium isotope fractionation during basalt differentiation: New measurements by multicollector sector ICP-MS, *Geochimica et Cosmochimica Acta* 63 907-910.

In addition to these offices, Thermo Electron Corporation maintains a network of representative organizations throughout the world.

Australia
+61 2 9898 1244

Austria
+43 1 333 50340

Belgium
+32 2 482 30 30

Canada
+1 800 532 4752

China
+86 10 5850 3588

France
+33 1 60 92 48 00

Germany
+49 6103 4080

Italy
+39 02 950 591

Japan
+81 45 453 9100

Netherlands
+31 76 587 98 88

Nordic
+46 8 556 468 00

South Africa
+27 11 570 1840

Spain
+34 91 657 4930

Switzerland
+41 61 48784 00

UK
+44 1442 233555

USA
+1 800 532 4752

www.thermo.com



Thermo Electron (Bremen) GmbH is certified DIN EN ISO 9001:2000

©2004 Thermo Electron Corporation. All rights reserved. All trademarks are the property of Thermo Electron Corporation and its subsidiaries.

Specifications, terms and pricing are subject to change. Not all products are available in all countries. Please consult your local sales representative for details.

AN30014_E 05/04C

Thermo
ELECTRON CORPORATION